



## Operating Instructions

### 1 Product Description

Use the Sequazyme™ BSA Test Standard Kit to test instrument function, optimize instrument parameters, and calibrate the mass scale using a known standard. The kit is designed for use in the Sequazyme™ Peptide Mass Standards Kit with the Voyager™ Biospectrometry™ Workstation or with any MALDI-TOF (matrix-assisted laser desorption ionization time-of-flight) mass spectrometry system.

The standard in the kit requires minimal preparation and covers a mass range of 60,000 to 70,000 daltons (Da). The standard was chosen for this kit on the basis of high purity and consistent results when analyzed using MALDI-TOF mass spectrometry.

#### Applications

Using the Sequazyme BSA Test Standard Kit, you can monitor results in a mass range from 60,000 to 70,000 Da for mass assignment, resolution, and sensitivity.

### 2 Materials

#### Materials Provided

The Sequazyme BSA Test Standard Kit includes —lyophilized bovine serum albumin, BSA (1 vial, 2 mg/vial)

#### Matrix and Diluent

- **Matrix B:Sinapinic Acid**—3,5-Dimethoxy-4-hydroxycinnamic acid, 1 vial (7 to 10 mg/vial)
- **Matrix B Diluent**—30% acetonitrile in 0.3% TFA, 1 vial (1 ml/vial)

After you use up the sinapinic acid matrix provided in the Sequazyme™ Peptide Mass Standards Kit, you can use commercially available sinapinic acid to analyze standards.

### 3 Preparing Matrix and Standards

See Section 7, Storing the Kit, for storage and stability conditions of prepared reagents.

**WARNING** CHEMICAL HAZARD. Diluent (with acetonitrile) is a flammable liquid and vapor. It may cause eye, skin, and respiratory tract irritation, central nervous system depression, and heart, liver, and kidney damage. Please read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

**CAUTION** CHEMICAL HAZARD. Sinapinic acid may cause eye, skin, and respiratory tract irritation. Please read the MSDS, and follow the handling instructions. Wear appropriate protective eyewear, clothing, and gloves.

#### Reconstitute BSA

Store the BSA Test Standard at –20° C prior to use. Sinapinic Acid Matrix from a Sequazyme kit is required to use this standard.

1. Add 0.5 ml HPLC-grade water to the BSA vial.
2. Recap and vortex for 1 minute.

You can use the standard immediately.

#### Sinapinic Acid Matrix

Use sinapinic acid matrix B from the Sequazyme Peptide Mass Standards Kit.

1. Add the entire contents of the Matrix B Diluent vial to the Matrix B:Sinapinic Acid vial.

Section	Page
1 Product Description.....	1
2 Materials.....	1
3 Preparing Matrix and Standards.....	1
4 Loading Matrix:Standard on Sample Plates.....	2
5 Analyzing Standards.....	2
6 BSA Standard Spectrum and Masses.....	2
7 Storing the Kit.....	2
8 Accessories, Spare Parts, and Ordering Information.....	3
9 Technical Support.....	3

- Vortex for 1 minute.
- Warm the solution, if necessary, to completely dissolve the solid.

Use the supernatant for sample preparation.

#### Standard

Final concentration of the standard after mixing with matrix is 4 pmol/μl.

- Pipette 1 μl of the reconstituted BSA solution into a microcentrifuge tube.
- Add 29 μl of the sinapinic acid matrix.
- Vortex briefly at low speed.

#### 4 Loading Matrix:Standard on Sample Plates

Each matrix:standard preparation yields 30 μl, enough to load approximately 30 sample positions.

- Pipette 1.0 μl of sample on the MALDI plate.
- Allow the mixture to air dry until all solvent is evaporated, usually less than 5 minutes.

Analyze within one day for best results.

#### 5 Analyzing Standards

##### Standard methods on your system

Your Voyager™ workstation includes standard methods created during factory testing of your instrument. Standard methods contain typical parameters for a given mass or mass range. Standard methods are useful as starting points when you create methods to analyze the standards in the Peptide Mass Standards Kit.

##### Analyzing standards

Analyze the standard according to the needs of your application.

Refer to Section 6, BSA Standard Spectrum and Masses, for representative spectra and mass assignments for standards and matrices in the kit.

#### 6 BSA Standard Spectrum and Masses

Figure 1 shows a representative spectrum for the test mixture. Masses are included in the spectra for peak identification only. Use the precise masses listed in Table 1 for calibration.

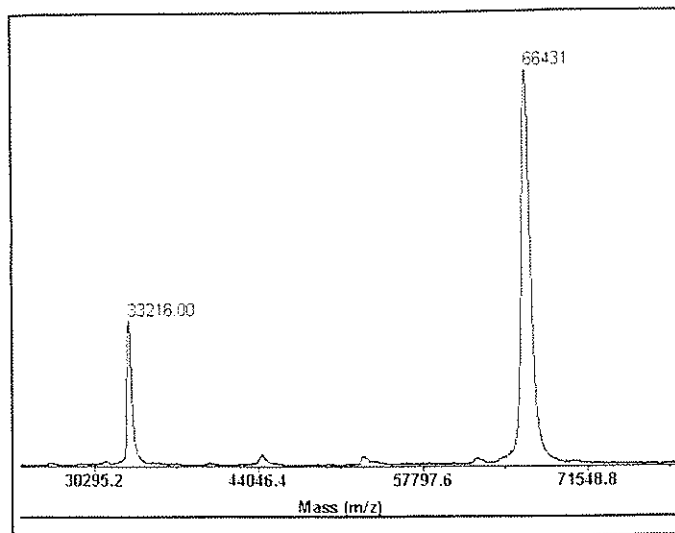


Figure 1 BSA Linear Resolution

Table 1 Mass Assignments for the Standard

Standard	Charge (n)	(M+nH) <sup>+</sup> Average
BSA (Bovine Serum Albumin)	+1	66,431
	+2	33,216

#### 7 Storing the Kit

Store the Sequazyme™ BSA Test Standard Kit and components of the kit under the following conditions. Avoid prolonged exposure to light.

Kit Component	Storage Temperature	Stability
Unopened kit	-20°C	1 year from date of shipment
Reconstituted standards	-20°C	6 months
Reconstituted matrix	4°C	1 week

## 8 Accessories, Spare Parts, and Ordering Information

Description	Quantity (Kits)	Part Number
Sequazyme™ Peptide Mass Standards Install Kit includes: <ul style="list-style-type: none"><li>• Sequazyme Peptide Mass Standards Kit</li><li>• BSA Test Standard Kit</li><li>• IgG1 Mass Standard Kit</li></ul>	1	4316866
Sequazyme Peptide Mass Standards Kit	1	P2-3143-00
BSA Test Standard Kit	1	2-2158-00
IgG1 Mass Standard Kit	1	GEN602151

## 9 Technical Support

PerSeptive Biosystems is committed to meeting the needs of your research through enabling technologies like the Sequazyme BSA Test Standard Kit. Our dedicated support staff is available to answer questions about using this product to the fullest extent possible.

PerSeptive Biosystems offers a complete line of time-of-flight (TOF) mass spectrometry products to meet your analysis needs. Please contact your PerSeptive Biosystems representative for technical and ordering information.


PerSeptive Biosystems publishes a continuing series of Application Notes and Technical Notes, highlighting ladder sequencing applications and strategies. Please contact PerSeptive Biosystems directly for a publication list.

For further details or for answers to questions related to other products contact PerSeptive Biosystems. Refer to the bottom of the following page for the number of the appropriate office.

Products and procedures described in this protocol are intended for research purposes only. Not for use in diagnostic procedures.

PerSeptive Biosystems and the fractal icon are registered trademarks and Sequazyme, Voyager, Biospectrometry, and Delayed Extraction are trademarks of PE Corporation or its subsidiaries in the U.S. and certain other countries.

Part Number 4317219, Rev. A, 06/00 ©2000 Printed In USA

 Printed on recycled paper.

---

PE Biosystems  
Headquarters  
850 Lincoln Center Drive  
Foster City, CA 94404  
Tel: (650) 638-5800  
Tel: (800) 345-5224  
[www.pebiosystems.com](http://www.pebiosystems.com)

PerSeptive Biosystems  
Division  
500 Old Connecticut Path  
Framingham, MA 01701  
Tel: (508) 383-7700  
Tel: (800) 899-5858  
[www.pbio.com](http://www.pbio.com)

For international office locations, please call the headquarters or refer to our web site at [www.pebiosystems.com](http://www.pebiosystems.com).

